

Cost-efficient single-use technology

3M™ Polisher ST is the next generation in single-use technology, designed as a cost-efficient option to increase biomass production in therapeutic recombinant protein processing. It is designed for use in most operating conditions for replacing the downstream Anion Exchange (AEX) polishing column in fed batch and/or continuous manufacturing processes. 3M Polisher ST is offered in a range of products suitable for early screening and process development of monoclonal antibody (mAb) processes and larger single use capsules for deployment in manufacturing processes in most polishing process conditions.

3M Polisher ST is a compact, small footprint solution. It has greater capacity for delivering higher purity and yield in the downstream polishing unit operation. This is a solution for a connected continuous downstream operation for the flexible facility of the future.

3M Polisher ST allows a facility strategy that can overcome scale limitations and enable cost-efficient manufacturing to support the growing demand for biologics.



3M Polisher ST is a synthetic, fully encapsulated single-use anion exchange device containing two complementary AEX-functional media: a quaternary ammonium ("Q") functional nonwoven (FNW) and a guanidinium (Gu)-functional membrane.

The Q-functional nonwoven provides reduction of turbidity (when present), DNA, and endotoxin, as well as a portion of the product's AEX capacity for host cell protein (HCP) and virus reduction.


The novel guanidinium functionality of the downstream AEX polishing membrane mimics the amino acid, arginine, one of the three positively charged naturally occurring standard amino acids that make up proteins. Comprising a resonance-delocalized positive charge and multiple co-planar pairs of hydrogen-bonding groups, the guanidinium group of arginine is observed in nature to interact with residues of the two negatively charged, carboxylic-acid functionalized amino acids, aspartic acid and glutamic acid, forming robust salt bridges that involve two hydrogen bonding interactions in addition to an electrostatic interaction.




Competitor

Small footprint of 3M Polisher ST (left) allows for more efficient use of space

3M Polisher ST

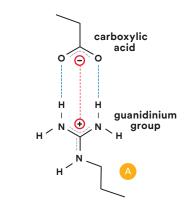

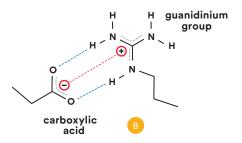


AEX performance of the guanidinium-functional membrane of 3M Polisher ST provides HCP and virus reduction in a salt tolerant (ST) manner compared with conventional Q-functional AEX resins and single-use media, which rely solely on electrostatic interactions that can be screened at elevated ionic strength.

Innovative design

The composite purification media comprises three primary components: Q-functional FNW, guanidinium-functional AEX membrane, and support membrane. The Q-functional portion of the AEX media is composed of four layers of polypropylene nonwoven that have been surface functionalized with a covalently attached quaternary ammonium functional polymer.

The guanidinium-functional portion of the AEX media is composed of three layers of a polyamide microporous membrane, having a 0.8-µm nominal pore size, that have been functionalized with a covalently attached guanidinium-functional polymer. The AEX media are followed by a single layer of a nominally 0.8-µm pore size polyamide membrane support.

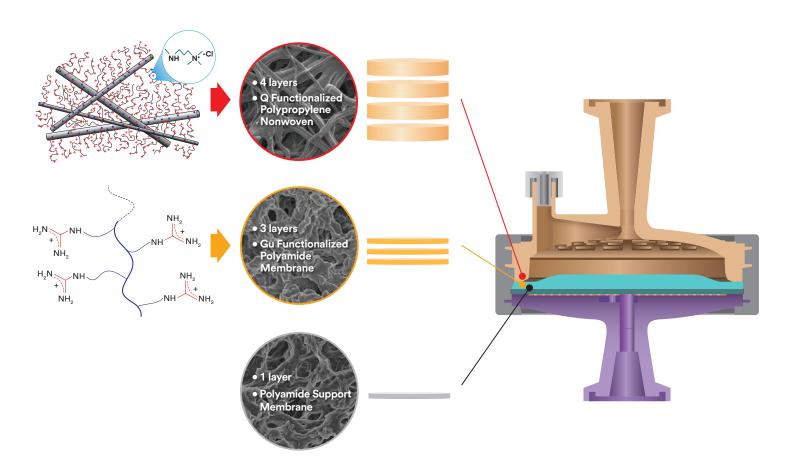

Functionalized membranes are used for high productivity chromatographic purification processes that match upstream efficiency without the need for oversizing.

3M Polisher ST single-use membrane adsorber purification solution is designed for the downstream processing of monoclonal antibodies (mAbs) and other recombinant proteins in flow-through mode.

3M Polisher ST is designed to improve mAb development and manufacturing processes in most polishing process conditions.

Purification – streamlined

Biopharmaceutical purification process improvements

This process train illustrates the potential of combining 3M products to work together to create an intensified manufacturing process, eliminating other process steps. 3M™ Zeta Plus™ **Depth Filter** • Removes whole cells and reduces cell debris • Sizes available from benchtop to production scale Master cell bank (MCB) Centrifuge 5000 Liter 50 Liter 500 Liter **Bioreactor** Bioreactor **Bioreactor** 3M™ Polisher ST • Fully encapsulated single-use AEX product Helps reduce HCP impurities, viruses and other negatively charged contaminants Intended for flow-through polishing chromatography Low pH Hold (Virus Inactivation) Column 3M™ LifeASSURE™ PDA Sterile Filter Polishing Column • 0.1 µm sterilizing membrane 3M™ Emphaze™ AEX Hybrid Purifier • Bacteria and mycoplasma removal • Robust turbidity reduction • DNA, HCP and endotoxin reduction 0.2 µm ЗМ™ Sterilizing Filter Filter R R R R

Performance

3M™ Polisher ST is designed for replacement of the Q-functional bead-based chromatography column with high throughput, cleaner effluent and lower total cost of ownership.

The nonwoven and membrane skeletons provide mechanical strength and durability, while the grafted hydrogel functionality creates a large three-dimensional surface area that contains a high density of functional groups with interconnected pores allowing for convective flow channels to achieve high flow rates during purification.

The high density of binding sites, together with a macroporous structure, enables high binding capacity for not only HCPs, but also large molecules, such as viruses and DNA. The high density of binding sites and low residence time results in high mAb recovery (>95%).

The high ligand density on our advanced anion exchange membrane provides robust impurity removal, viral clearance, and a 50 to 100X higher mAb loading capacity than chromatography resin beads to further reduce media volume requirement. The impurity clearance performance

(HCP ≤1000 ppm, ≤100 ppb DNA, and viral clearance) is independent of the load which could be as high as 10 kg/m², enabling downsizing of the unit operation.

3M Polisher ST offers a combination of high dynamic antibody-loading (10 kg/m²) capacities with low residence time (about 0.2 min compared to 1 minute for traditional Q resin), opening the way to an intensified high-productivity, truly single-use purification platform.

Robust viral clearance is achieved over a range of buffer and pH conditions, which provide a wide design space for operation, as shown in the data below. Performance is maintained at high conductivity (20 mS/cm) even in citrate buffer, which reduces the need for feed dilution or buffer exchange requirement resulting in a simplified process with reduced cost factors.

AEX chromatography removes HCP and DNA and is also an effective virus removal step as part of an orthogonal viral clearance technology platform. Under process conditions of pH 5.5 to 7.5 and 5 to 20 mS/cm, 4 to 7 LRV can be achieved using 3M Polisher ST.




Test Data

Dynamic binding capacity


3M™ Polisher ST lab and scale-up capsules have target mAb loading capacities from 1 g to 1 kg.

Larger capsules with target mAb loadings of 2.25 kg and 16 kg are planned. Scalability of capsules with different frontal media surface area was assessed by measurement of bovine serum albumin (BSA) protein dynamic binding capacity (DBC) as a simulant for HCP. The BSA DBC per unit area is quite constant across all capsules, having a mean value of 22.2 g/cm² and a standard deviation of 0.98 g/cm².

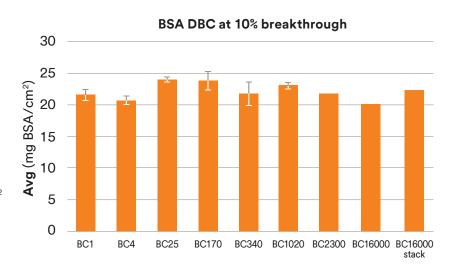


Consistent DBC under a wide range of pH and conductivity

The BSA DBC specification for the functional membrane only is 7.8-12 mg BSA/cm² in 25 mM Tris, pH 8, 50 mM NaCl. High salt (20 mS/cm) reduces the capacity of the membrane by ~2-3 mg/cm². Additional capacity provided by the FNW at conductivities < ~10 mS/cm.

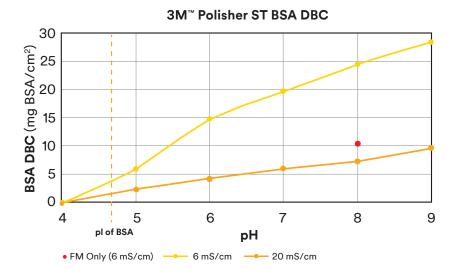


The differential pressure ($\triangle P$) specification for the functional membrane only is < 10 psid in 25 mM Tris, pH 8, 50 mM NaCl at 1 mL/min/cm². Lower salt (< 6 mS/cm) slightly increases pressure due to contribution from FNW. Maximum differential pressure (all capsule sizes) is 35 psid.

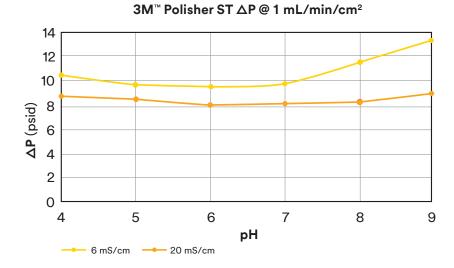

HCP reduction


- ~500 ppm HCP in a VIN mAb pool
- > 60% removal in Phosphate buffer,
- > 22% removal in Citrate buffer

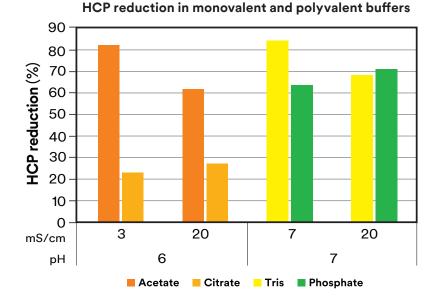


In comparison to Q column


3M™ Polisher ST offers a combination of high dynamic antibody-loading (10 kg/m²) capacities with low residence time (about 0.2 min compared to 1 minute for Q-resin), opening the way to an intensified high-productivity, truly single-use purification platform.

Chromatographic columns have a limited loading capacity for polishing and virus capture. Furthermore, they are large and have a high investment cost. Traditional column chromatography systems are slow in producing recombinant proteins.

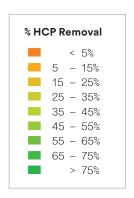

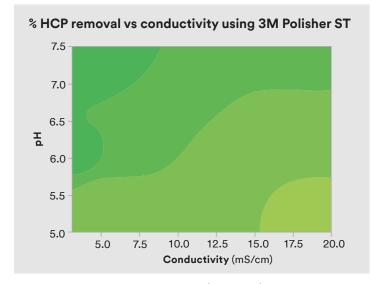



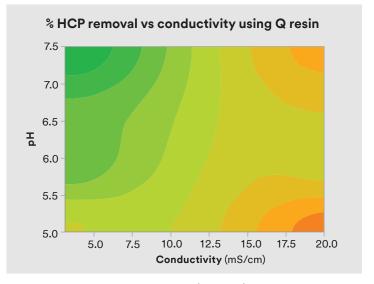



Q Chemistry (200 g/L)


Test data (continued)

mAb recovery

Without significant adjustments to pH and conductivity, 3M™ Polisher ST has robust performance with respect to HCP and DNA clearance and better mAb recovery.

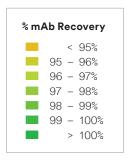

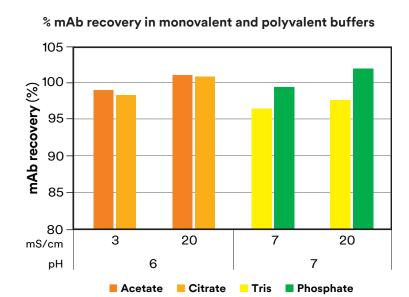


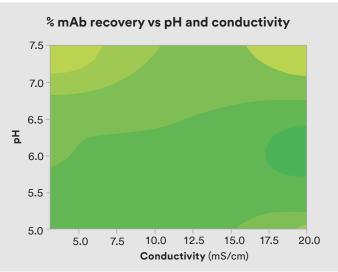

3M Polisher ST (10 kg/m²) in acetate/tris

DNA reduction

DNA clearance across a variety of initial concentrations (~10-1000 ppb) down to detection limit levels (0.05 ppb DNA) in all buffers (acetate, citrate, phosphate, tris) and conductivities were tested and are reported in the table alongside.

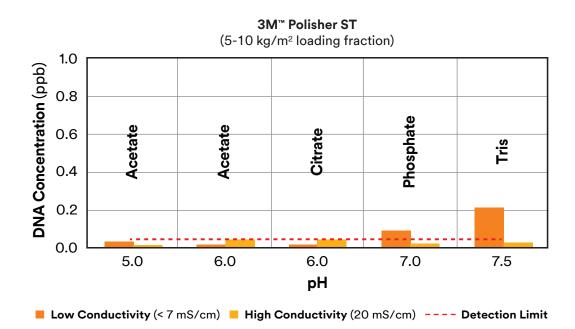

HCP and DNA reduction and mAb recovery in legacy, modern and next gen mAb feed streams

Buffer	рН	Conductivity (mS/cm)	Feed stream	Turbidity (NTU)	Target HCP (ppm)	Target DNA (ppb)	HCP reduction (%)	mAb recovery (%)	DNA conc. (ppb)	DNA reduction (LRV)
Tris	7.5	5	Legacy	no	200	20	81.1	102	0.05	2.7
Phosphate	7.5	20	Legacy	no	200	20	58.2	100	< 0.05	> 2.9
Acetate	5.5	5	Modern	yes	500	1000	60.8	100	< 0.05	> 5.8
Citrate	5.5	5	Modern	no	500	50	18.5	99.6	0.06	3.5
Tris	7.0	7	Next-gen	no	500	50	81.7	99.7	< 0.05	> 3.1
Acetate	6.0	7	Next-gen	no	500	50	75.0	100	< 0.05	> 3.7

Load 10 kg/m²

Performance in turbid streams

3M Polisher ST was challenged with turbid streams to simulate different downstream processes.

Buffer	рН	Conductivity (mS/cm)	Turbidity (NTU)	Initial mAb conc. (mg/mL)	Initial HCP conc. (ppm)	mAb recovery (%)	HCP reduction (%)	Fraction	DNA conc. (ppb)	DNA reduction (LRV)
Acetate-Tris	5.5	6	39.7	22.3	1465	98.6	80	Load	47560	
								0-5 kg/m²	< 0.05	> 5.98
								5-10 kg/m ²	0.09	5.73
Tris-Acetate	7.0	6	18.3	23.4	1508	100.2	54	Load	106800	
								0-5 kg/m ²	2.18	4.69
								5-10 kg/m ²	3.34	4.50

Similar HCP (> 50%) and DNA reduction (> 4 LRV) and mAb recovery (> 97%) compared to non-turbid samples. The ΔP across the capsule did not significantly change during the 10 kg/m² load.

Legacy

Post Depth Filter and CEX Bind and Elute step. Depth Filter and CEX have reduced some contaminant levels.

Post VIN with minimal pH adjustment. Process in general is cleaner compared to legacy but there is less preconditioning prior to AEX resulting in higher contaminant and turbidity levels.

Next-gen

Similar feed stream to modern with less DNA and turbidity and slightly higher pH and conductivities for subsequent polishing steps (HIC/MM).

Test data (continued)


Viral Clearance in Buffer


Viruses XMuLV, Reo-3, PrV and MVM


XMuLV, Reo-3 and PrV all showed > 6 LRV clearance (detection limit) from pH 5 - 7.5, conductivities 3 - 20 mS/cm and in both monovalent (Acetate/Tris) and polyvalent (Citrate/Phosphate) buffers.


MVM showed > 4 LRV clearance from pH 5.5 - 7.5, conductivities 5 - 20 mS/cm and in both monovalent (Acetate/Tris) and polyvalent (Citrate/Phosphate) buffers.

Viral Clearance studies were performed with MVM and XMuLV using mAb solutions representing legacy, modern and next-gen mAb feed streams. The Tris pH 7.5, 5 mS/cm MVM study was performed at Charles River. All other studies were performed at Texcell.

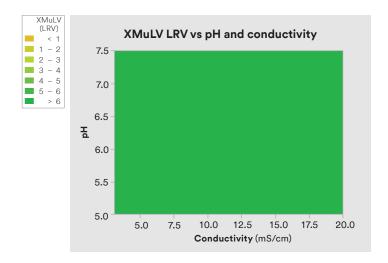

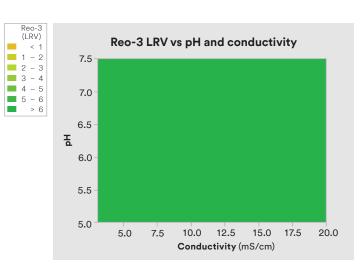

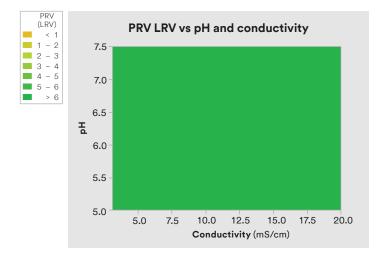

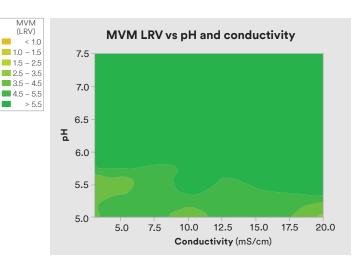



Buffer	рН	Conductivity (mS/cm)	Feed stream	Turbidity (NTU)	Target HCP	Target DNA	Avg viral clearance (LRV)			
							MVM		XMuLV	
					(ppm)	(ppb)	0-5 kg/m²	5-10 kg/m ²	0-5 kg/m²	5-10 kg/m ²
Tris	7.5	5	Legacy	no	200	20	6.16	6.53	> 6.26	> 6.26
Phosphate	7.5	20	Legacy	no	200	20	> 4.69	> 7.41	> 6.61	> 6.89
Acetate	5.5	5	Modern	yes	500	1000	5.34	5.86	> 6.53	> 7.54
Citrate	5.5	5	Modern	no	500	50	5.33	6.14	6.16	4.74
Tris	7.0	7	Next-gen	no	500	50	5.60	6.05	> 6.88	6.41
Acetate	6.0	7	Next-gen	no	500	50	> 5.43	5.92	> 6.72	> 7.72

3M™ Polisher ST is designed as a next generation bioprocessing solution to address the challenge biopharmaceutical companies have with the adoption of new manufacturing methods, for continued growth and innovation. Within the next five years, it is estimated that approximately 35 percent of today's biologics will be manufactured using some sort of process intensification methods. Facilities will be smaller but capable of producing higher volumes of multiple molecules.

The key to this advancement will be the use of methods such as continuous bioprocessing technologies, some of which may be single use, which would enable facilities to reduce both their size and cycle times. The complexity of the bioprocessing industry has challenged companies' innovation efforts, due to potential costs required for changes to take place.

3M Polisher ST is a single-use technology designed as a cost-efficient option to increase biomass production in therapeutic recombinant protein processes in all operating conditions and to replace downstream AEX polishing column in fed batch and/or continuous manufacturing.

3M Polisher ST is a compact, small footprint, higher capacity solution for delivering higher purity and yield in the downstream polishing unit operation in a connected continuous downstream operation for the flexible facility of the future.

3M Polisher ST enables a facility strategy that can overcome scale limitations and enable cost-efficient manufacturing to support the growing demand for biologics.





Intended Use: 3M[™] Polisher ST single-use filter products are intended for use in biopharmaceutical processing applications of aqueous based pharmaceuticals (drugs) and vaccines in accordance with the product instructions and specifications, and cGMP requirements, where applicable. Since there are many factors that can affect a product's use, the customer and user remain responsible for determining whether the 3M product is suitable and appropriate for the user's specific application, including user conducting an appropriate risk assessment and evaluating the 3M product in user's application.

Prohibited Use: 3M advises against the use of these 3M products in any application other than the stated intended use(s), since other applications have not been evaluated by 3M and may result in an unsafe or unintended condition. Do not use in any manner whereby the 3M product, or any leachable from the 3M product, may become part of or remains in a medical device that is regulated by any agency, and/or globally exemplary agencies, including but not limited to: a) FDA, b) European Medical Device Directive (MDD), c) Japan Pharmaceuticals and Medical Devices Agency (PMDA) or in applications involving permanent implantation into the body; Life-sustaining medical applications; Applications requiring food contact compliance.

Product Selection and Use: Many factors beyond 3M's control and uniquely within user's knowledge and control can affect the use and performance of a 3M product in a particular application. As a result, end-user is solely responsible for evaluating the product and determining whether it is appropriate and suitable for end-user's application, including a completing a risk assessment that considers the product leachable characteristics and its impact on drug safety, conducting a workplace hazard assessment and reviewing all applicable regulations and standards (e.g., OSHA, ANSI, etc.) Failure to properly evaluate, select, and use a 3M product and appropriate safety products, or to meet all applicable safety regulations, may result in injury, sickness, death, and/or harm to property.

Warranty, Limited Remedy, and Disclaimer: Unless a different warranty is specifically stated on the applicable 3M product packaging or product literature (in which case such warranty governs), 3M warrants that each 3M product meets the applicable 3M product specification at the time 3M ships the product. 3M MAKES NO OTHER WARRANTIES OR CONDITIONS, EXPRESS OR IMPLIED, INCLUDING, BUT NOT LIMITED TO, ANY IMPLIED WARRANTY OR CONDITION OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR ARISING OUT OF A COURSE OF DEALING, CUSTOM, OR USAGE OF TRADE. If a 3M product does not conform to this warranty, then the sole and exclusive remedy is, at 3M's option, replacement of the 3M product or refund of the purchase price.

Limitation of Liability: Except for the limited remedy stated above, and except to the extent prohibited by law, 3M will not be liable for any loss or damage arising from or related to the 3M product, whether direct, indirect, special, incidental, or consequential (including, but not limited to, lost profits or business opportunity), regardless of the legal or equitable theory asserted, including, but not limited to, warranty, contract, negligence, or strict liability.



3M Purification Inc.
3M Separation and Purification Sciences Division

400 Research Parkway Meriden, CT 06450 USA

Phone: 1-800-243-6894 1-203-237-5541 Web: 3M.com/polisher 3M, Emphaze, LifeASSURE and Zeta Plus are trademarks of 3M Company. All other trademarks are property of their respective owners.

© 2021 3M Company. All rights reserved. Please recycle. Printed in U.S.A. 70-2016-0021-3